

SNP Analysis 1: DNA Extraction

Serhat Sevli, C. Yunus Sahan

Abstract



Nehir Biyoteknoloji Ltd. www.nehirbt.com

Citation: Serhat Sevli, C. Yunus Sahan: SNP Analysis 1: DNA Extraction. **protocols.io**

10.17504/protocols.io.gycbxsw

Published: 21 Jan 2017

Before start

This protocol is derived from

[Reference: Genematrix Quick Blood DNA Purification Kit, version 1.2, EURx company]

Protocol

Activation of column

1. Apply 40µl of activation Buffer QB onto the spin-column (do not spin) and keep it at room temperature till transferring lysate to the spin-column.

*Warning #1

Addition of Buffer QB onto the center of the resin enables complete wetting of membranes and maximal binding of DNA.

*Warning #2

The membrane activation should be done before starting isolation procedure.

Sample preparation

2. Add 200µl blood/body fluid into a new 1.5-2 mL tube.

Add 10µl Proteinase K and next 200µl Sol QB buffer.

Vortex the mixture thoroughly and incubate for 20 min at 60°C inside an incubator.

Add 200µl 96% ethanol in the same tube of DNA sample.

Vortex the mixture thoroughly and centrifuge for 1 min at 12000 rpm.

*Warning #3

For sample volumes less than 200µl, add PBS to adjust the volume to 200µl.

*Warning #4

If RNA-free DNA is crucial for downstream applications, add 2µl RNase A. Mix by vortexing and incubate 5 min at room temperature.

*Warning #5

If purifying DNA viruses, it is recommended to start with 200µl serum or plasma to prepare pure viral DNA (cellular DNA-free)

Washing of lysate DNA

3. Transfer the lysate sample into the spin-column and place in a new clean collection tube.

Centrifuge for 2 min at 12000 rpm.

Take out the spin-column, discard flow-through and place back the spin-column in the collection tube.

Add 500µl Wash QBX1 buffer to the spin-column and centrifuge for 1 min at 12000 rpm.

Take out the spin-column, discard flow-through and place back the spin-column in the collection tube.

Add 500µl Wash QBX2 buffer to the spin-column and centrifuge for 2 min at 12000 rpm.

* **Warning #6**

Continue centrifugation at all steps, if not all of the sample passed through the column.

Collection of DNA

4. Place the spin-column in a new collection tube (1.5-2 mL)

Add 50-200 µl of Elution buffer (10 mM Tris-HCl, pH 8.5) heated to 60°C to elute bound DNA in column.

Incubate the spin-column/collection tube assembly for 5 min at room temperature with the cap is closed.

Centrifuge the spin-column for 1 min at 12000 rpm.

* **Warning #7**

Addition of the elution buffer directly onto the center of the resin improves DNA yield. To avoid transferring traces of DNA between the spin-column do not touch the spin column walls with the micropipette.

* **Warning #8**

The researcher is free to use other types of Elution solutions regarding further use of DNA sample.

* **Warning #9**

The elution can be repeated once again as described in steps 15-17. This step improves DNA recovery from the column. A new collection tube can be used to prevent dilution of the first eluate or collection tube from step 15 can be reused to combine the eluates.

* **Warning #10**

More than 200µl should not be used to elute into a single 1.5 ml microcentrifuge tube, as the spin-column will come into contact with the eluate, causing DNA contamination.

Storage of DNA sample

5. Discard the spin-column, cap the collection tube, label as appropriate.

DNA is ready for further experiment

It can be stored either at 2÷8°C or at -20°C.

Warnings

- * All the related steps must be done in a biotech lab using appropriate clothing and equipment.
- * All centrifuging steps must be performed by an expert.
- * Please learn well about the hazards of all chemical used in this protocol.